



# Acquisition Protocol for Cytek® cFluor® MDSC Kit on PBMCs

## Contents

Introduction .....	1
Preparing SpectroFlo® Software .....	2
Add fluorochromes to the library .....	2
Import the experiment template.....	2
Setting up the Instrument.....	3
Acquiring Controls and Samples .....	5
Acquire controls in Reference Group .....	5
Unmix reference controls .....	5
Acquire multicolor samples .....	7
Appendix A: Reusing Single Color Controls.....	10
Appendix B: Single Color Control Signatures for 3-Laser (V-B-R) Cytek® Northern Lights™ Suggested Gating When Using Cells as Controls.....	11
Appendix C: Example of Similarity™ Matrix for Cytek® cFluor® MDSC Kit .....	14
Appendix D: Example of Adjusted Spillover Matrix .....	15
Appendix E: Example of Gating in Multicolor Sample on PBMCs .....	16

## Introduction

This acquisition protocol provides step-by-step instructions to set up your Cytek® Northern Lights™ system (3-laser V-B-R configuration or higher) for data acquisition of the Cytek® cFluor® MDSC Kit. This protocol provides instructions on 1) preparing SpectroFlo® software, 2) setting up the instrument, and 3) acquiring controls and samples.

This kit contains 25 tests.

**NOTE:** For suggestions on how to prepare human peripheral mononuclear cells and whole blood, see **Sample Preparation (PBMCs) Protocol for Cytek® cFluor® MDSC Kit**

## Preparing SpectroFlo® Software

### Add fluorochromes to the library

1. Import **Cytek® MDSC Kit Tags\_PBMC**.csv file into the **Library** module, under **Fluorescent Tags** to add the fluorochromes used in this kit.

Name	Created By	Date Created	Date Modified	Emission Wavelength	Laser Excitation	Display Name	Synonyms
Blue Laser	System	May 12, 2021 - 12:36 PM	March 02, 2023	cFluor B515	515	488	cFluor B515
Red Laser	System	May 12, 2021 - 12:36 PM	March 02, 2023	cFluor B548	548	488	cFluor B548
Violet Laser	System	May 12, 2021 - 12:36 PM	March 02, 2023	cFluor G575	575	488	cFluor G575
Fluorescent Proteins	System	May 12, 2021 - 12:36 PM	May 12, 2021	cFluor BYG610	610	488	cFluor BYG610
Viability	System	May 12, 2021 - 12:36 PM	May 12, 2021	cFluor BYG667	667	488	cFluor BYG667
cFluor Immunoprofiling Kit Tags, 25 Color	Admin	June 24, 2021 - 10:42 AM	June 24, 2021	cFluor BYG710	710	488	cFluor BYG710
MDSC Kit Tags_Blood	Admin	March 17, 2023 - 11:09 AM	March 17, 2023	cFluor BYG781	781	488	cFluor BYG781
MDSC Kit Tags_PBMC	Admin	March 17, 2023 - 11:09 AM	March 17, 2023	cFluor R659	659	635	cFluor R659
				cFluor R685	685	635	cFluor R685
				cFluor R720	720	635	cFluor R720
				cFluor R840	840	635	cFluor R840
				cFluor V450	450	405	cFluor V450
				cFluor V505	505	405	cFluor V505
				Ghost Dye V540	540	405	Ghost Dye V540

**NOTE:** If any fluorochromes were previously entered into SpectroFlo® Library, a warning message will appear. Click **OK** to overwrite the information in the library with the new fluorescent tag information.


### Import the experiment template

1. **Cytek® MDSC Kit Template\_PBMC** includes a reference group with predefined stopping criteria, assigned marker names, as well as recommended acquisition and analysis worksheets.
2. Import the template into the **Library** module, under **Experiment Templates**.


Name	Created By	Date Created	Description	Shared
Cytek MDSC Kit Template_PBMC	Admin	March 14, 2023 - 10:04 AM		<input type="checkbox"/>
Cytek MDSC Kit Template_Blood	Admin	March 14, 2023 - 10:04 AM		<input type="checkbox"/>
Default	Admin	March 02, 2023 - 16:47 PM	Default Experiment	<input type="checkbox"/>
14C IP Panel QC Template	Admin	November 19, 2021 - 10:21 AM		<input type="checkbox"/>
Cytek 8C TBMNK QC Template	Admin	November 19, 2021 - 10:15 AM		<input type="checkbox"/>
cFluor Immunoprofiling Kit, 25 Color Template	Admin	June 24, 2021 - 10:47 AM	Default Experiment	<input type="checkbox"/>

## Setting up the Instrument

1. Follow the instructions for instrument setup and **Performing Daily QC** as outlined in the User's Guide.
2. From the **Acquisition** module, create a new experiment by clicking on **Template**, then choose "**Cytek® MDSC Kit Template\_PBMC.**" Click **Open** and give a name to your experiment.

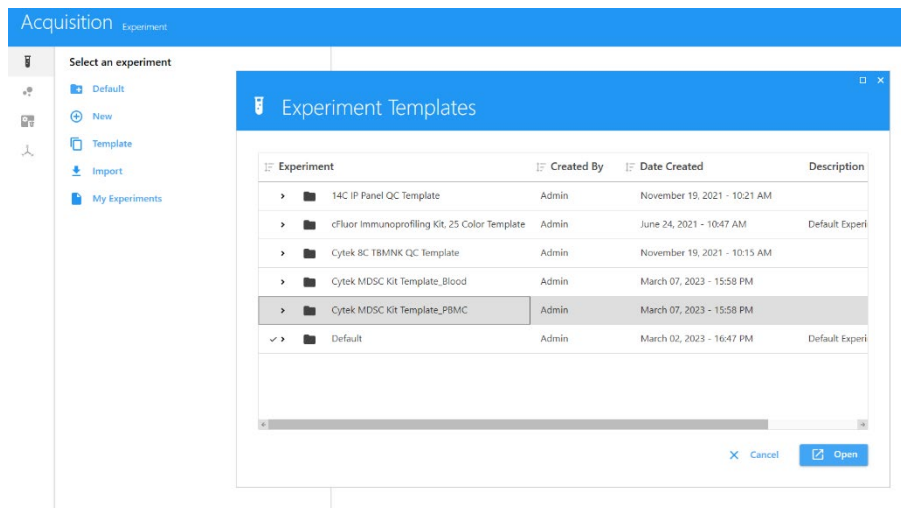

3. For reference controls it is recommended to run cells for single color controls unless otherwise stated. **NOTE:** See Table 1 for reference control type recommendations for each marker.
4. Each sample tube is set to acquire certain number of cells. This can be changed by clicking **Edit**, then changing the Stopping Gate and Stopping Criteria under **Acquisition**.

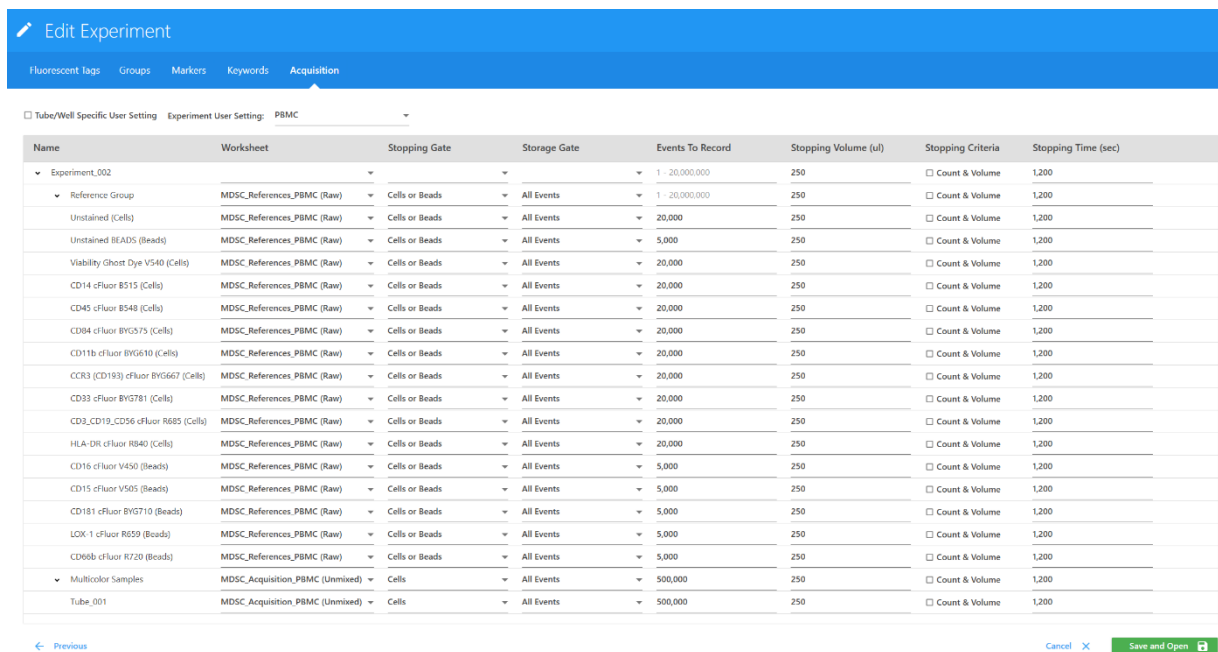

5. Add tubes and groups as needed. To preserve the predefined acquisition conditions, duplicate the existing tubes or groups.

**Table 1.** Reference Control Type Recommendations

<b>Laser</b>	<b>Target</b>	<b>Fluorochrome</b>	<b>Recommended Control Type</b>
Violet	CD16	cFluor® V450	Beads
	CD15	cFluor® V505	Beads
	Viability	Ghost Dye™ V540	Cells
Blue	CD14	cFluor® B515	Cells or Beads
	CD45	cFluor® B548	Cells or Beads
	CD84	cFluor® BYG575	Cells or Beads
	CD11b	cFluor® BYG610	Cells or Beads
	CCR3 (CD193)	cFluor® BYG667	Cells or Beads
	CD181	cFluor® BYG710	Beads
	CD33	cFluor® BYG781	Cells or Beads
Red	LOX-1	cFluor® R659	Beads
	CD3	cFluor® R685	Cells or Beads
	CD19		
	CD56		
	CD66b	cFluor® R720	Beads
	HLA-DR	cFluor® R840	Cells or Beads


**NOTE:** Recommendations are for use with Cytex® FSP™ CompBeads.

## Acquiring Controls and Samples

### Acquire controls in Reference Group


1. Preview unstained cell control at low flow rate to minimize wasted sample volume. Starting from the default CytekAssaySetting, optimize the FSC and SSC gains, as well as the threshold to fully visualize the cells of interest (see Figure 1).

**Note:** Instrument settings can be saved for future use by clicking the **Save As** button in the Instrument Control window. The gains for all fluorescent parameters are set up with Cytek Assay Setting in the instrument and only FCS, SSC and threshold need to be optimized for specific sample types.

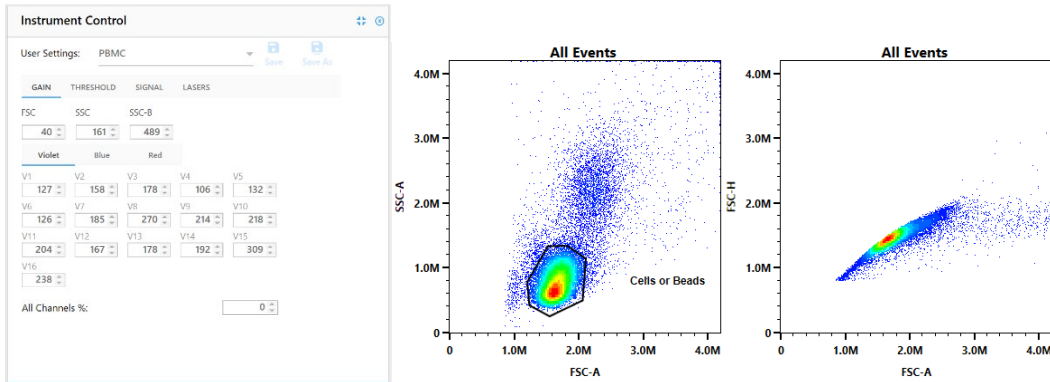

**Figure 1:** Example cell resolution of human PBMC samples after FSC and SSC adjustment with threshold set at FSC 50,000.

2. Acquire unstained and single color controls in the Reference Control group using the **MDSC\_References\_PBMC** worksheet. Click **Start** to preview for 5 to 10 seconds until the event rate stabilizes, then click **Record** to record each sample.

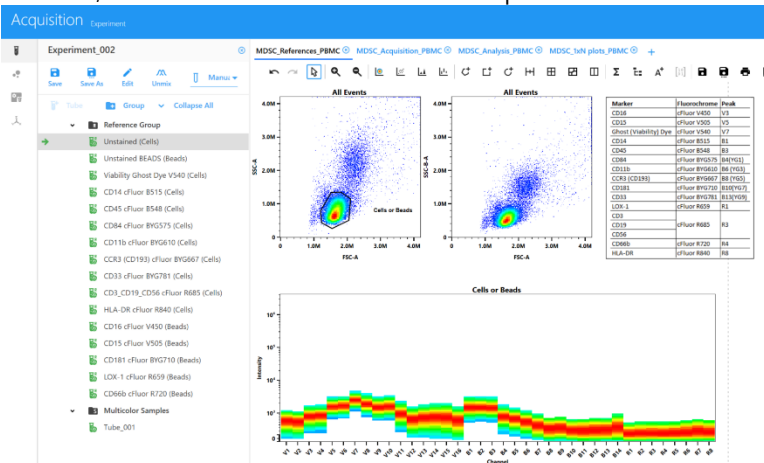


### Unmix reference controls

1. Once all controls have been acquired, Click **Unmix**.  
**NOTE:** Refer to Appendix A for additional workflows to store and reuse the reference controls.
2. Under **Select Controls** tab in the Unmix Experiment wizard, ensure that:
  - 1) Under **UNSTAINED CONTROLS** "Use Control from Experiment" is selected, and the additional **Unstained BEADS** control is included.

- 2) Unstained is selected as **Reference Group – Unstained (Cells)** for all markers stained on cells except for the following which should use marker internal negative signal as negative reference: Ghost Dye V540, CD84, CD3/CD19/CD56, HLA-DR, and CCR3 (CD193).
- 3) For markers LOX-1, CD16, CD15, CD66b, and CD181 Unstained is selected as **Reference Group – Unstained BEADS (Beads)**
- 4) **Autofluorescence as a Fluorescence Tag** is selected for 5 laser instruments and unselected for 3 laser instruments. Click **Next**.

**Note:** Autofluorescence extraction is ALWAYS recommended for data generated in 5-laser systems. For data generated in 3-laser systems (V-B-R), there is a need to assess the level of autofluorescence of the sample by checking the spectrum plot of unstained cells. If autofluorescence is low in unstained cells, as shown in Appendix B, autofluorescence extraction is NOT recommended.

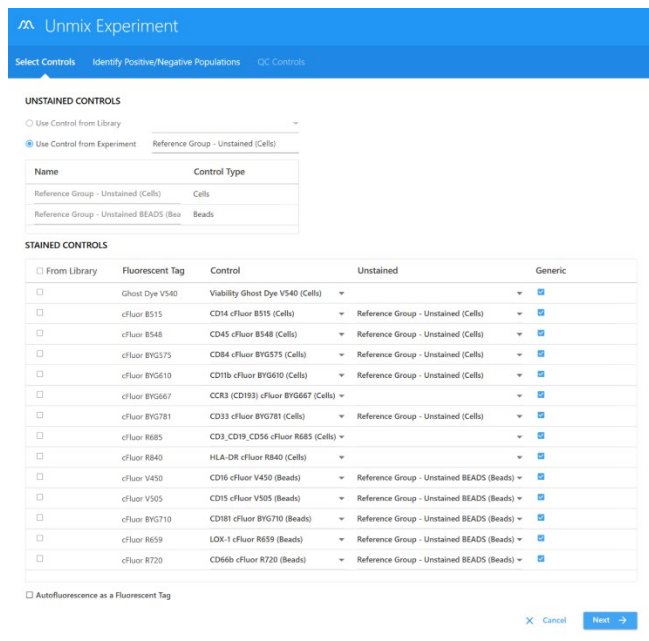

**UNSTAINED CONTROLS**

Use Control from Library

Use Control from Experiment: Reference Group - Unstained (Cells)

Name	Control Type
Reference Group - Unstained (Cells)	Cells
Reference Group - Unstained BEADS (Beads)	Beads

**STAINED CONTROLS**


<input type="checkbox"/> From Library	Fluorescent Tag	Control	Unstained	Generic
<input type="checkbox"/>	Ghost Dye V540	Viability Ghost Dye V540 (Cells)		<input type="checkbox"/>
<input type="checkbox"/>	cFluor B515	CD14 cFluor B515 (Cells)	Reference Group - Unstained (Cells)	<input type="checkbox"/>
<input type="checkbox"/>	cFluor B548	CD45 cFluor B548 (Cells)	Reference Group - Unstained (Cells)	<input type="checkbox"/>
<input type="checkbox"/>	cFluor BYG575	CD84 cFluor BYG575 (Cells)	Reference Group - Unstained (Cells)	<input type="checkbox"/>
<input type="checkbox"/>	cFluor BYG610	CD11b cFluor BYG610 (Cells)	Reference Group - Unstained (Cells)	<input type="checkbox"/>
<input type="checkbox"/>	cFluor BYG667	CCR3 (CD193) cFluor BYG667 (Cells)		<input type="checkbox"/>
<input type="checkbox"/>	cFluor BYG781	CD33 cFluor BYG781 (Cells)	Reference Group - Unstained (Cells)	<input type="checkbox"/>
<input type="checkbox"/>	cFluor R685	CD3, CD19, CD56 cFluor R685 (Cells)		<input type="checkbox"/>
<input type="checkbox"/>	cFluor R840	HLA-DR cFluor R840 (Cells)		<input type="checkbox"/>
<input type="checkbox"/>	cFluor V450	CD16 cFluor V450 (Beads)	Reference Group - Unstained BEADS (Beads)	<input type="checkbox"/>
<input type="checkbox"/>	cFluor V505	CD15 cFluor V505 (Beads)	Reference Group - Unstained BEADS (Beads)	<input type="checkbox"/>
<input type="checkbox"/>	cFluor BYG710	CD181 cFluor BYG710 (Beads)	Reference Group - Unstained BEADS (Beads)	<input type="checkbox"/>
<input type="checkbox"/>	cFluor R659	LOX-1 cFluor R659 (Beads)	Reference Group - Unstained BEADS (Beads)	<input type="checkbox"/>
<input type="checkbox"/>	cFluor R720	CD66b cFluor R720 (Beads)	Reference Group - Unstained BEADS (Beads)	<input type="checkbox"/>

Autofluorescence as a Fluorescent Tag

3. Under **Identify Positive/Negative Populations** tab, ensure for all single color controls that:
  - 1) Scatter plot is gated on the appropriate population,
  - 2) Black bar is on the peak channel,
  - 3) Signature of each fluorochrome matches the expected spectrum, and
  - 4) Negative and positive gates in the histogram are correctly positioned.

**Note:** Refer to Appendix B for the correct gate positioning, expected spectra and peak channels of each fluorochrome, and the positioning of negative and positive gates in the histograms.

3. Click **Next**. Under the **QC Controls** tab, click on **Similarity Matrix** to confirm all controls were appropriately stained. Click on **View Similarity Index** to compare the expected complexity index value found in Appendix C.

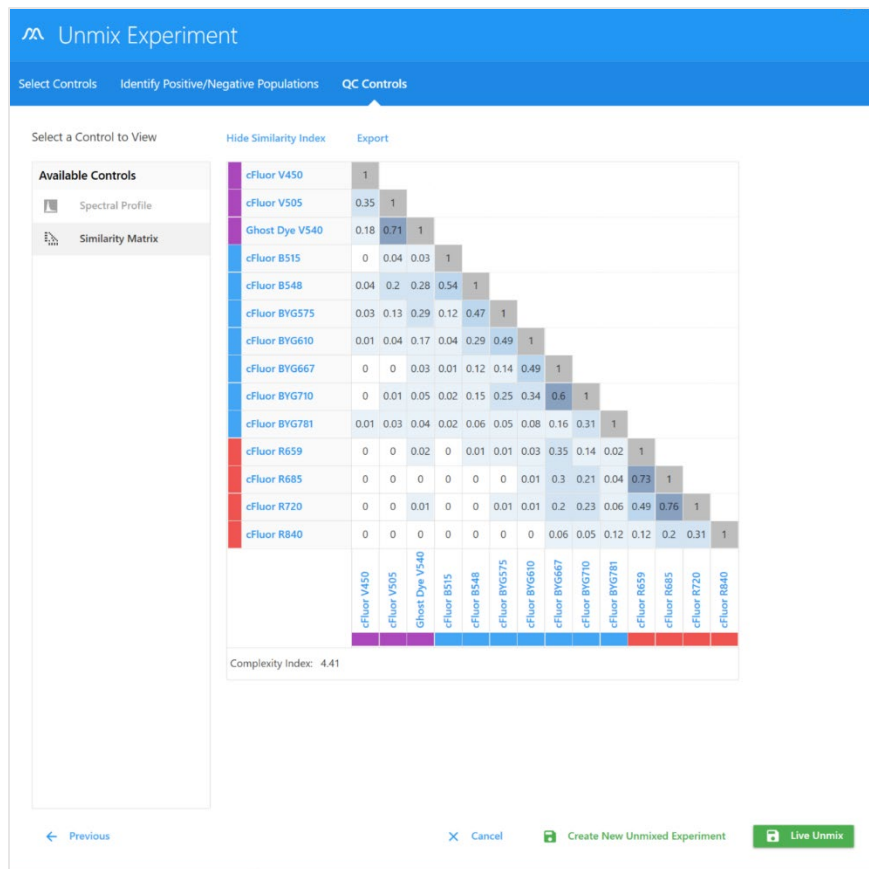

4. Click **Live Unmix**.

### Acquire multicolor samples


1. To acquire multicolor samples, **MDSC\_Acquisition\_PBMC** worksheet should be selected in the Acquisition tab to collect multicolor samples in the experiment template.
2. Ensure the same settings for the FSC and SSC gains used in the Unstained control are used for the multicolor samples. These FSC and SSC gains can be verified by right clicking on the tube in the experiment, selecting **Edit Properties** and checking on the **Instrument Settings** tab. If needed, the scatter gains can be adjusted.

### Analyze multicolor samples

3. A **MDSC\_Analysis\_PBMC** worksheet is included in the experiment template to be used for multicolor sample data analysis. In this worksheet, an additional Time Plot is included as an initial plot to eliminate any abnormal cells (i.e. abnormal cells due to edge effect).
4. Also included in the experiment template is a **MDSC\_1xN\_plots\_PBMC** worksheet. Select and acquired multicolor sample and open this worksheet to view the multicolor sample in the worksheet.
5. In case the recorded multicolor samples need spillover adjustment, right click on the multicolor tube. From the drop-down menu, click **Edit Properties**.





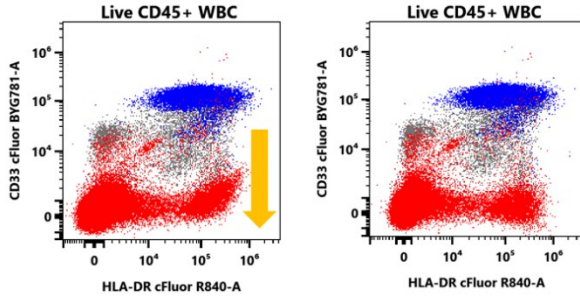

**Figure 2:** Example of unmixing error adjustment.


9. Click **Save** and close out of the Tube/Well Properties wizard. The adjusted compensation matrix can be applied to other multicolor samples by copying and pasting Spillover, if similar unmix errors are observed.
10. Repeat the unmixing error adjustment for all multicolor samples as needed
11. Manually adjusted compensation matrix can be saved to library for future use if using SpectroFlo® version 3.0 or higher.

**NOTE:** Refer to an example of adjusted spillover matrix in Appendix D.

12. Double click on the multicolor to open **MDSC\_Analysis\_PBMC** worksheet to view and adjust gates in the worksheet.

**NOTE:** Refer to Appendix E for an example of a PBMC multicolor sample.

13. To export an experiment:
  - 1) Save and close the experiment,
  - 2) Click **My Experiments**,
  - 3) Select the experiment and click **Export**, and
  - 4) Choose a directory and click **Export**.

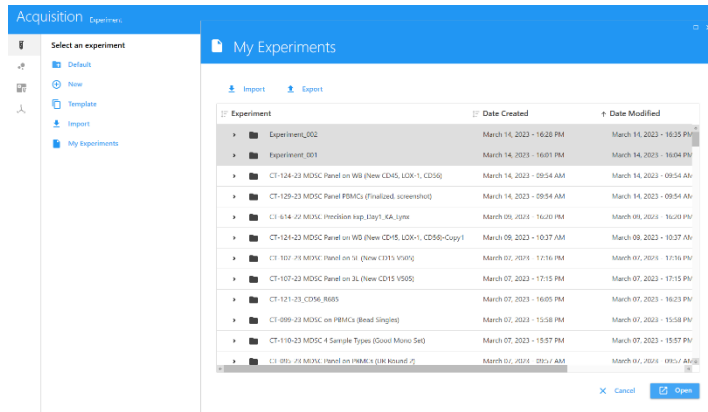


## Appendix A: Reusing Single Color Controls

To reuse reference controls in future experiments, follow the instructions below.

**NOTE:** For best results, the instrument should be maintained properly, QC performed daily and the same reagent lot used. The samples need to be collected in the same instrument as the reference controls.


### Reusing single color controls from a previous experiment


1. From the **Acquisition** module, open **My Experiments**. Right click the saved experiment and select **Duplicate**. This will duplicate the experiment with the reference controls and multicolor samples.
2. Under Multicolor Group, right click on the sample to duplicate or add tubes for the new multicolor samples
3. Delete any old multicolor samples that were carried over with data
4. Click **Edit** to open Edit Experiment wizard
5. Under **Groups**, right click on Multicolor and select **Unstained Control** to add a group specific unstained control for Multicolor Sample

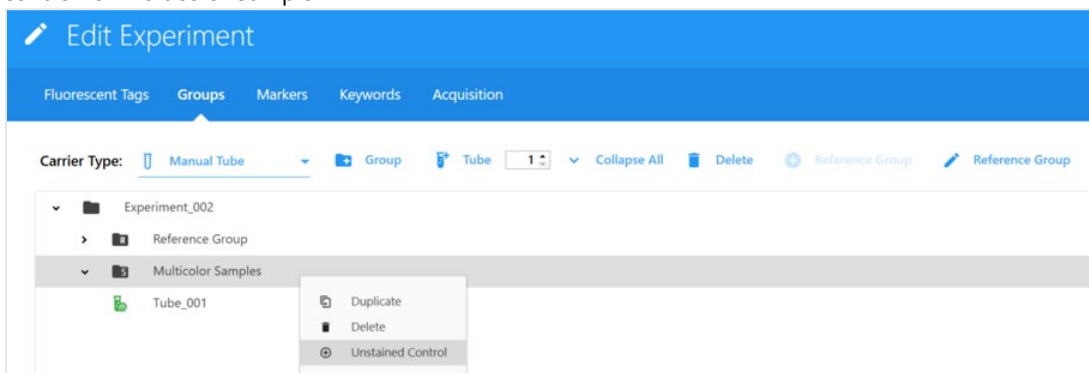


6. Under **Acquisition**, define Worksheet, Stopping Gate, and Events to Record
7. Click **Save and Open**
8. Preview and record the unstained control
9. Live Unmix with **Autofluorescence as a Fluorescence Tag** selected for 5 laser instruments and unselected for 3 laser instruments. Click **Next**.  
*Note: Autofluorescence extraction is ALWAYS recommended for data generated in 5-laser systems. For data generated in 3-laser systems (V-B-R), there is a need to assess the level of autofluorescence of the sample by checking the spectrum plot of unstained cells. If autofluorescence is low in unstained cells, as shown on page 13, autofluorescence extraction is NOT recommended.*
10. Acquire multicolor samples
11. Adjust the gates in the **MDSC\_Analysis\_PBMC** worksheet as needed.

## Appendix B: Single Color Control Signatures for 3-Laser (V-B-R) Cytex® Northern Lights™ Suggested Gating When Using Cells as Controls

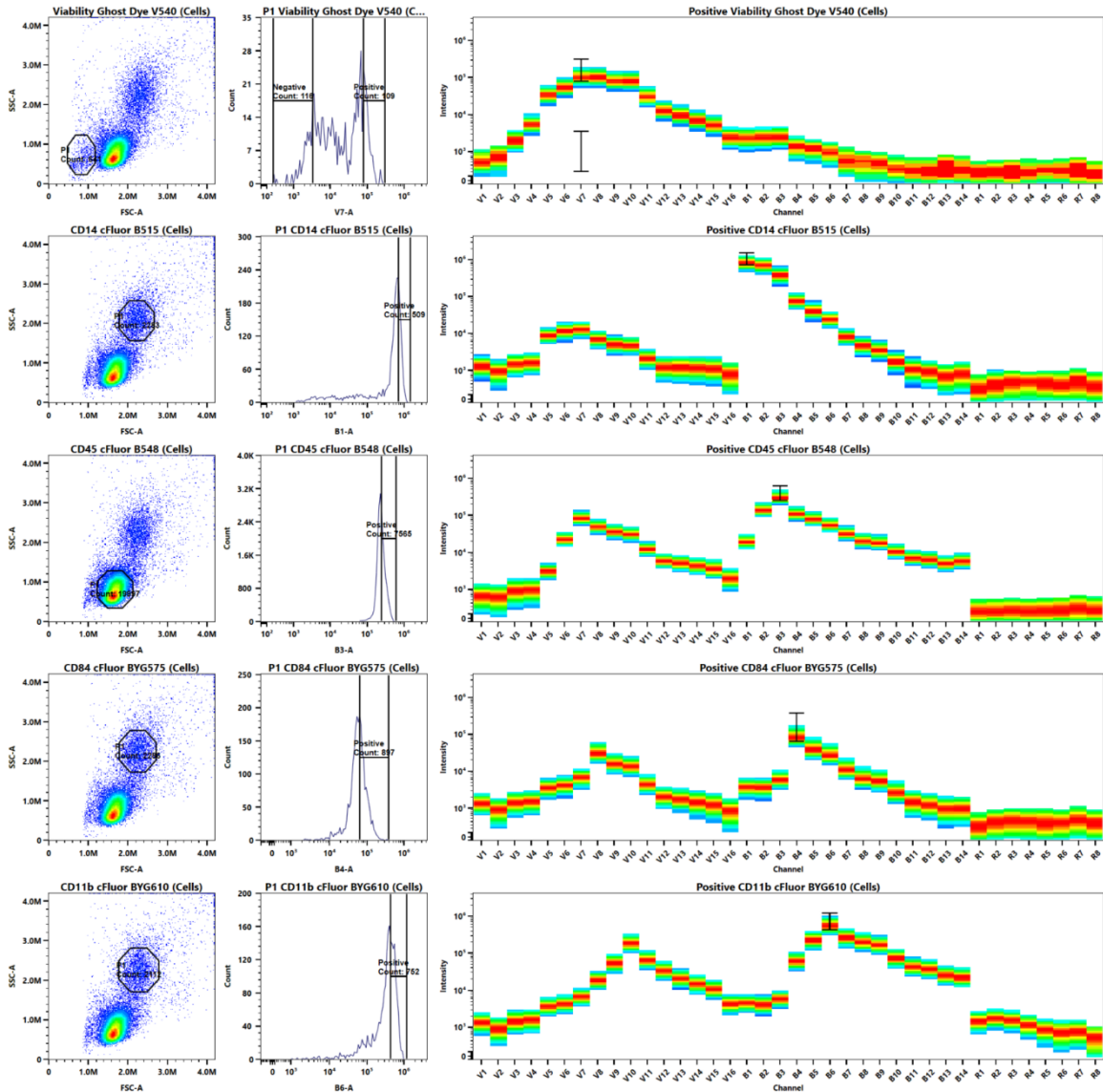

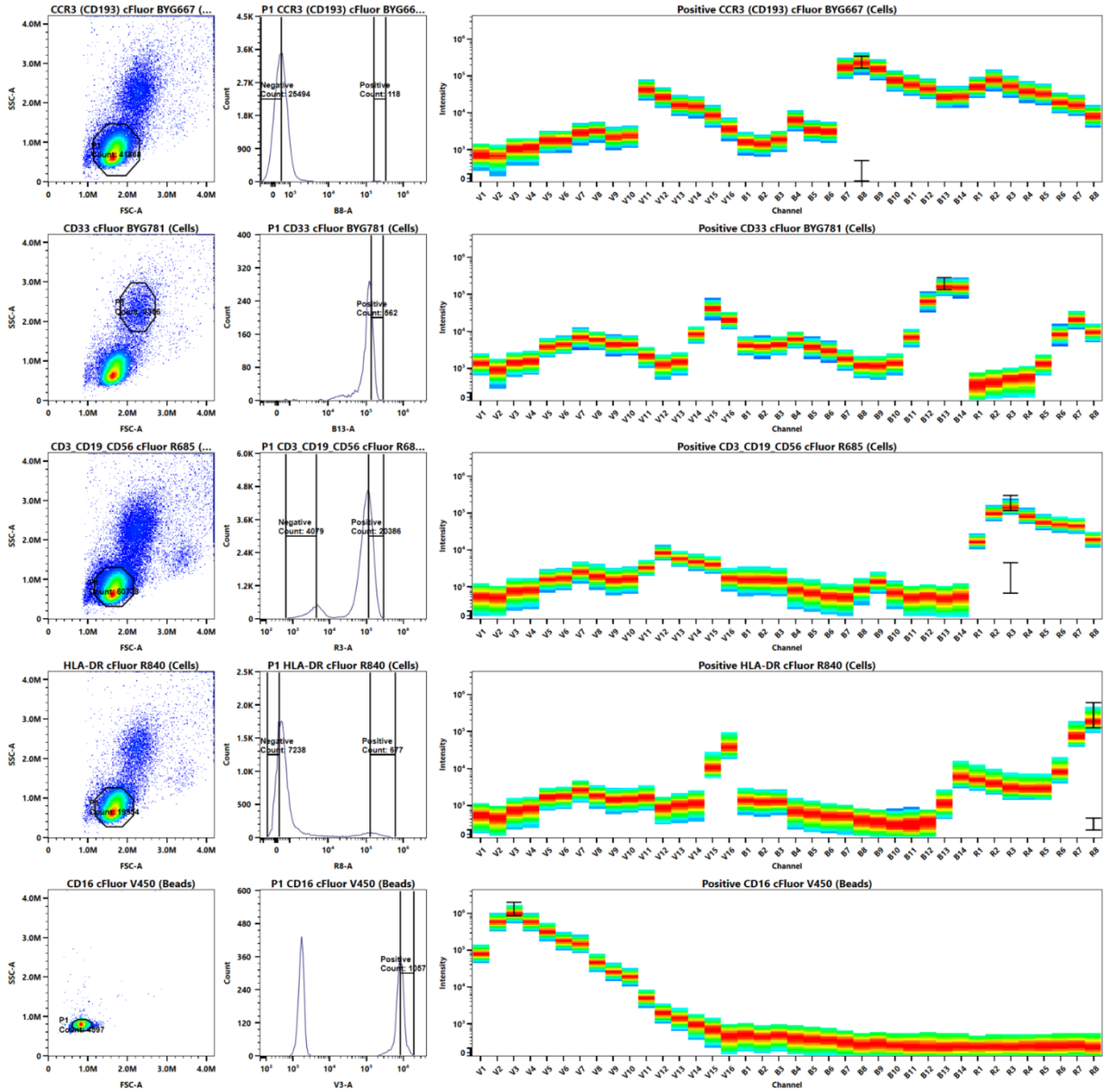

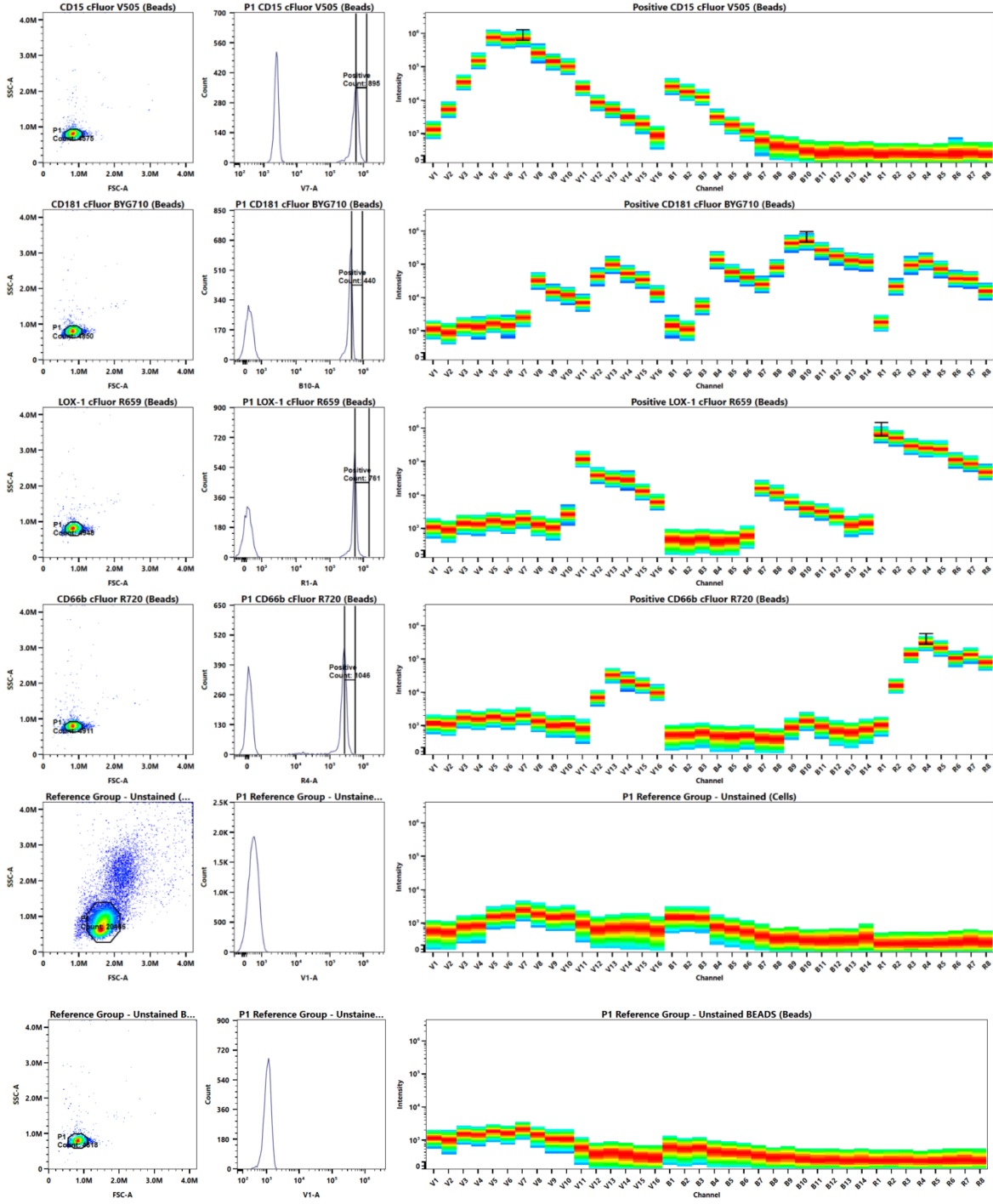



## Appendix D: Example of Adjusted Spillover Matrix

Tube/Well Properties

Tube/Well    Fluorescent Tag    Acquisition    Keywords    Instrument Settings    Spillover

Enable Compensation   
 Apply From Library   
 Apply From File   
 Save To Library   
 Save To File

INTO + FROM +	cFluor V450	cFluor V505	Ghost Dye V540	cFluor B515	cFluor B548	cFluor BYG575	cFluor BYG610	cFluor BYG667	cFluor BYG710	cFluor BYG781	cFluor R659	cFluor R685	cFluor R720	cFluor R840
cFluor V450	100	0	0	0	0	0	0	0	0	0	0	0	0	0
cFluor V505	0	100	0	-0.5	0	0	0	0	0	0	0	0	0	0
Ghost Dye V540	0	0	100	0	0	0	0	0	0	0	0	0	0	0
cFluor B515	0	0	0	100	0	0	0	0	0	0	0	0	0	0
cFluor B548	0	0	0	0	100	0	0	0	0	0	0	0	0	0
cFluor BYG575	0	0	0	0	0	100	0	0	0	0	0	0	0	0
cFluor BYG610	0	0	0	0	0	0	100	0	0	0	0	0	0	0
cFluor BYG667	0	0	0	0	0	0	0	100	0	0	0	0	0	0
cFluor BYG710	0	0	0	0	0	0	0	0	100	0	0	0	0	0
cFluor BYG781	0	0	0	0	0	0	0	0	0	100	0	0	0	0
cFluor R659	0	0	0	0	0	0	0	0	0	0	100	0	0	0
cFluor R685	0	0	0	0	0	0	0	0	0	0	0	100	0	0
cFluor R720	0	0	0	0	0	0	0	0	0	0	0	0	100	0
cFluor R840	0	0	0	0	0	0	0	0	0	0	0	0	0	100

**Figure 4:** Corrections made by checking NxN permutation plots for 14 fluorochromes. Please note that all needed corrections are below 3%. If corrections larger than 5% need to be applied, carefully QC your reference controls, comparing the staining pattern and gating positioning provided in Appendix B.







**For Research Use Only. Not intended for use in diagnostic procedures.**

cFluor® B515, cFluor® B548, cFluor® R685, cFluor® R720, and cFluor® R840 are equivalent to CF® 488A, CF® 514, CF® 660C, CF® 700, and APC-CF® 790T respectively, manufactured and provided by Biotium, Inc. under an Agreement between Biotium and Cytek (LICENSEE). The manufacture, use, sale, offer for sale, or import of the product is covered by one or more of the patents or pending applications owned or licensed by Biotium. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product for the purchaser's own internal research. No right under any other patent claim, no right to perform any patented method, and no right to perform commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is conveyed expressly, by implication, or by estoppel.

Cytek® FSP™ CompBeads are developed and manufactured by Slingshot Biosciences, Inc., Emeryville, CA.

"Cytek", "Northern Lights", "FSP", "cFluor", "Ghost Dye", "SpectroFlo", "Similarity" and "Complexity" are trademarks or registered trademarks of Cytek Biosciences, Inc. All other service marks, trademarks and tradenames appearing herein are the property of their respective owners.